




**Biomedical Marketing Associates**  
**Wexford, PA**  
**www.BiomedicalMarketing.com**  
**phone: (724) 935-6840**

**Retrieving Tool**  
**BCT-101**  
**\$95.00 ea.**

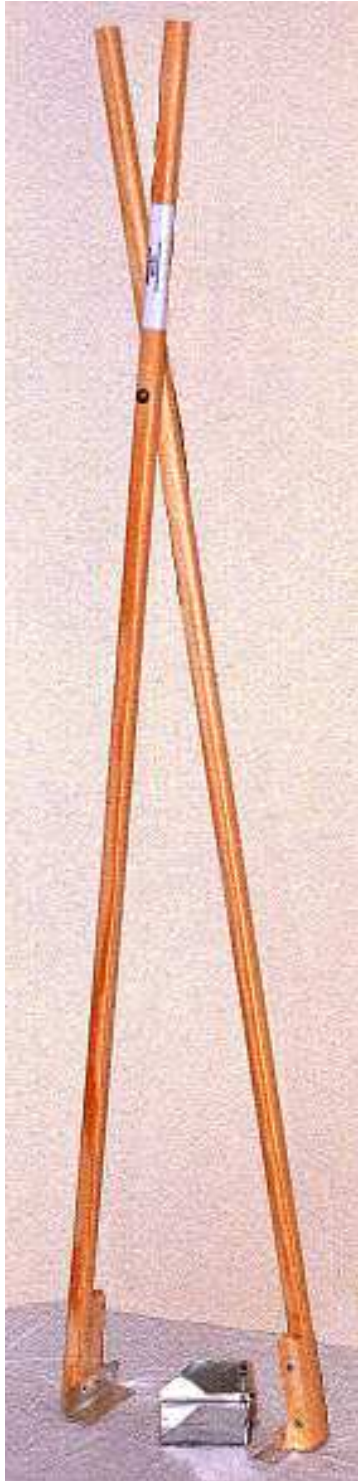

Designed to help recover objects dropped into liquid nitrogen freezers. What if you drop a sample tube, a pen, or even a whole storage box into a cryogenic storage vessel. How do you recover it?

At 4' long, the **Retrieving Tool** can be used in most liquid nitrogen freezers with neck openings greater than 5" for safe recovery of lost items. The wooden probes can be immersed in liquid nitrogen without chilling the handle area.

There is a detachable steel basket for recovering individual smaller items. For larger items, such as boxes, the basket is removed to leave 2 metal angle feet to support the item being recovered.

*Never* put your hand into liquid nitrogen. *Never* put your head below the rim of a large open top freezer. The dangers are too great for frostbite and asphyxiation. Use the **Retrieving Tool** and stay safe.



Send purchase orders to  
Biomedical Marketing Associates  
by fax (724) 935-6840  
by email: [sales@BiomedicalMarketing.com](mailto:sales@BiomedicalMarketing.com)

prices & specifications subject to change without notice  
freight prepaid and added to invoice